

## Sample Preparation for SampleJet

- Buy New 7” Tubes from Storeroom with Purple “Biospin” Caps
- 8” Tubes Will Not Fit in Carrousel
- Use 1 – 30 mg of Sample (10-30 for  $^{13}\text{C}$ )
- Minimum Solvent Depth 3.5 cm!
- Top of Tube Level and Smooth, No Cracks
- Cap Grips Tube Firmly, Not Loose
- No Labels of Any Kind On Tube: You May Write with Felt Pen Just Below Cap
- Enter Each Sample in Log to Keep Track