

Bruker “Open Shop” Automation Steps

- Insert 7” Tube With Special Cap in Rack: Holders 1-96 Around Periphery of Carrousel
- Write Login Name, Holder Number, Sample ID and Date in Log
- Login to ICON-NMR Automation Window
- Double-Click on Holder in Upper Half of Window
- Enter Sample ID (Filename), Expno., Experiment
- Add Any Additional Experiments
- Submit to Queue
- Optional: Check Progress of Queue at <http://aviii400.chem.arizona.edu:8015>
- When Finished, Pick Up Zip File of Data at <http://aviii400.chem.arizona.edu>
- Pick Up Your Sample From Carrousel or Overflow Racks